

Dialysis Tubing Preparation

1. Cut tubing into 18-24 inch lengths.
2. Curl into a flask solution of 2% sodium bicarbonate/ 1mM EDTA.
3. Boil for 10 minutes.
4. Rinse 2X with dH₂O.
5. Boil for 10 minutes in a solution of 1 mM EDTA.
6. Allow to cool and store at 4°C in a fresh solution of 1mM EDTA in sterile dH₂O.

SOLUTIONS

Sodium bicarbonate solution:

1000mls dH₂O
20 g sodium bicarbonate
2 mls EDTA