

Revised 3/2002
Julie McLear

Ethanol Precipitation of Probe

Elute probe as usual (quiagen kit)
Save 1 μ l for scintillation count

Start with 100 μ l probe (or other volume)
Add 10 μ l 3 M NaAc, 250 μ l EtOH, 1-2 μ l tRNA
Mix solution
Incubate at -20 C overnight
Spin 15 min. @ 4 C
Remove supernatant
Wash/add 500 μ l 70% EtOH
Spin 15 min
Remove supernatant
Air dry the pellet
Resuspend in 50 μ l 1 X TE pH 8.0 (or dH₂O)
Take 1 μ l for scintillation count
Compare probe before /after column and before/after precipitation